

## SLT Series Science Lab Table

### Leg Assembly Instructions

The SLT Series has been designed with the capability to be used with both our standard non-book compartment style or with book compartments (purchased separately).

To assemble the table using the standard non-book compartment style, follow the below instructions.

#### Tools Required

WRENCH



**Step 1:** Place the tabletop upside-down on the floor with apron side up.

*Note: Floor should be clear of any objects or dirt in order not to damage the tabletop.*

**Step 2:** Remove plastic covers on bolts and place each leg in position with leg bolts going through the corner braces.

**Step 3:** Place a washer, lock washer and nut onto each bolt and tighten well using the provided wrench.

**Step 4:** Slide a leg boot onto each leg assuring that the wider side is facing the bottom of the leg.

**Step 5:** Screw a glide into the bottom of each leg.

**Step 6:** Turn the table over with the tabletop facing up. Place the table in desired location. Adjust the glides to assure that table is stable.

**Step 7:** Slide down protecting boots to cover the adjusting glides.

*Note: Extra apron marked BC is for use with book compartments only.*

---

Dear Customer,

Thank you for your purchase. Although extreme care goes into manufacturing & packaging of this product, mistakes are possible. If you encounter any missing parts or difficulty in assembly. etc. Please call our customer service **TOLL FREE at 1-800-261-4112** for courteous & immediate resolution to your problem.

**CAUTION:** Double check all screws to be tightened securely before use. Check all screws periodically.